



For young plant culture test

Water-soluble TANEPIITA!

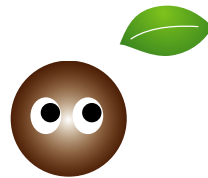


On the round water-soluble sheet, adhesive dots are placed in same interval (5 x 4). TANEPIITA saves your time and labor for seeding, and helps to execute the germination test in stable condition.

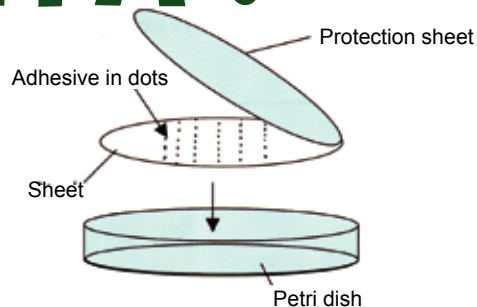

How to use

1. Peel off the protection sheet.
2. Place the sheet on petri dish, etc. and a pinch of seeds.
3. Spread the seeds over the adhesive, tilting the dish left and right.
4. After checking that 20pcs of seeds are stuck in order, hold TANEPIITA! with tweezers.

Example for young plant culture test

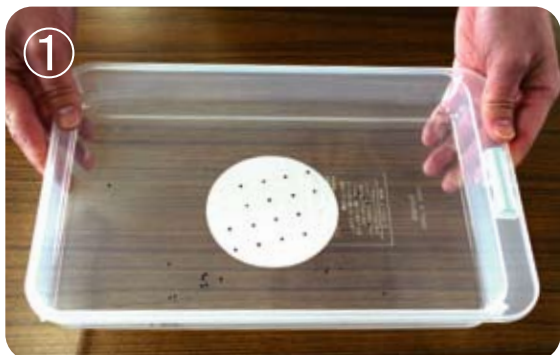

Peel off the protection sheet of water-soluble TANEPIITA! and sprinkle a pinch of seeds on adhesive.

※If necessary, correct the seeds in order manually.



Remove TANEPIITA! from the dish and place it on pot.



Cover the water-soluble TANEPIITA ! with 20g soil and power appropriate volume of soil.



Specifications

Sheet dia. : 90mm, water-soluble paper
Dot : 20/sheet, interval 16mm,
dia. of dot 1.0mm
Adhesive : Acrylic adhesive
Protection paper : Releasing paper
Recommended environment : Approx.23°C Humidity 65%
Validity : Recommended to use in 6months
Unit : 100pcs/box

Water-soluble

TANEPITA !



Cautions

Please read before starting to use.

【Cautions for use】

- Take care not to collect dust or touch by fingers on the adhesive dots. Otherwise the seeds will not be stucked.
- It is possible that more than 1 pc of seed is stucked on a dot. Please adjust by tweezers, etc.
- Do not handle by wet hands.
- When the room temperature is low, adhesive could be harden and loose some viscosity.
- Appropriate environment for use is temperature 23°C, humidity approx.65%.

【Cautions for storage】

- Recommended period for storage is up to 6 months.
- Keep out of direct sunlight and store at 3 ~ 4°C.
- Keep in plastic bag to avoid high humidity or drying.
- Protection paper shall not be peeled off from adhesive.

14-3-1000S

FHK Fujihira Industry Co., Ltd.

6-11-6 Hongo Bunkyo-ku Tokyo 113-0033 Japan

Tel: +81-3-3812-2270 Fax: +81-3-3812-3663

E-mail : overseas@fujihira.co.jp

URL: <http://www.fujihira.co.jp>